


**Instruction Sheet for
the PASCO
Model ME-9495A**


ANGLE INDICATOR

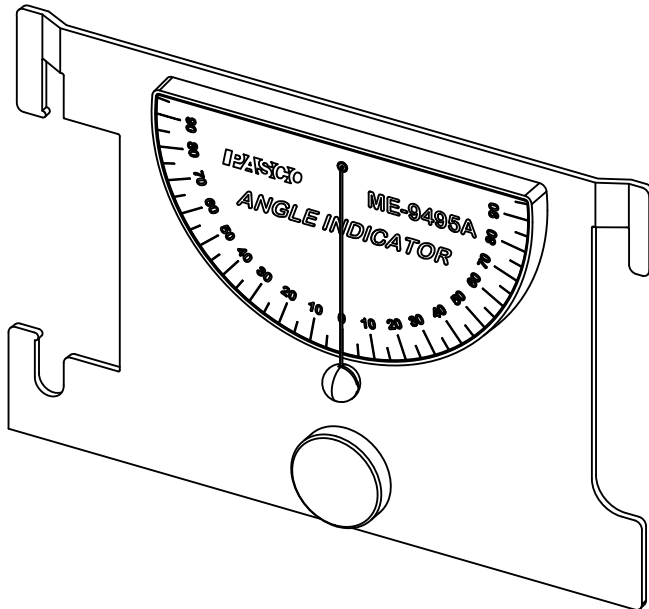


Introduction

The PASCO ME-9495A Angle Indicator is an accessory used for measuring angles from 0° to 90° in two directions. There are four different methods of using the Angle Indicator:


- hand-held, by suspending the device over a string.

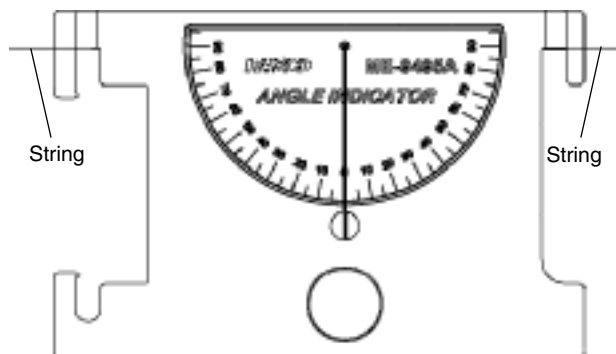

- attached to a PASCO Dynamics Cart, by sliding the device into the slots on the top of the cart.

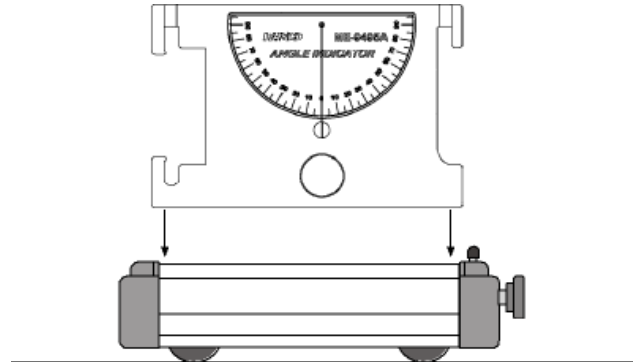


- attached to the side of a PASCO Dynamics Track.

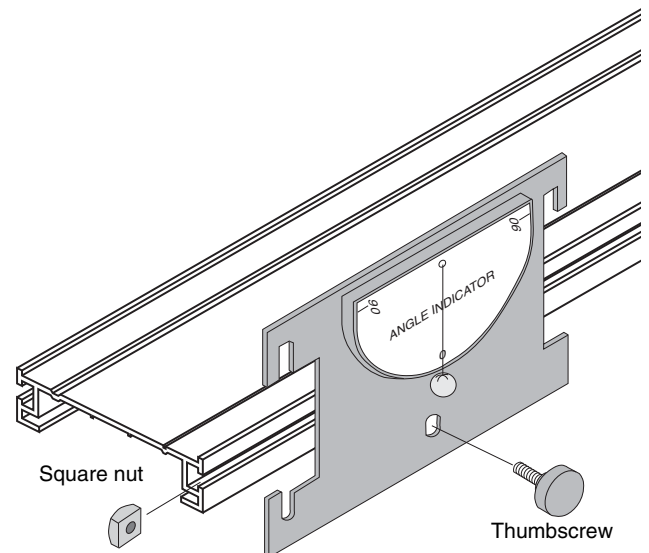


© 2009 PASCO scientific

- slipped over the platform on the PASCO Rotary Motion Sensor.

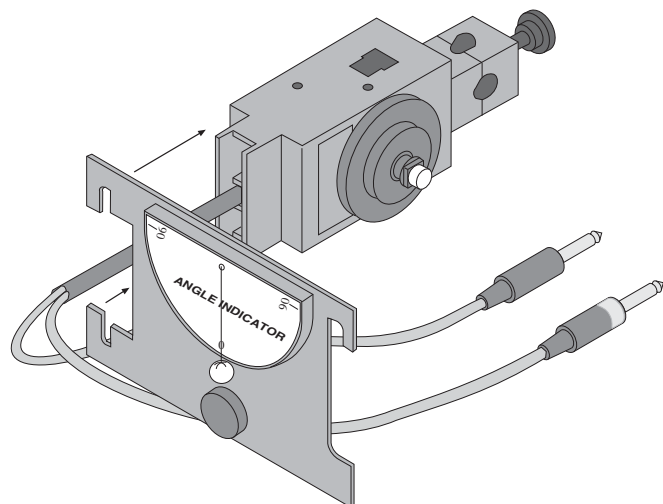

Maintenance

The thread that supports the spherical weight is attached to the back of the Angle Indicator with a hex screw.

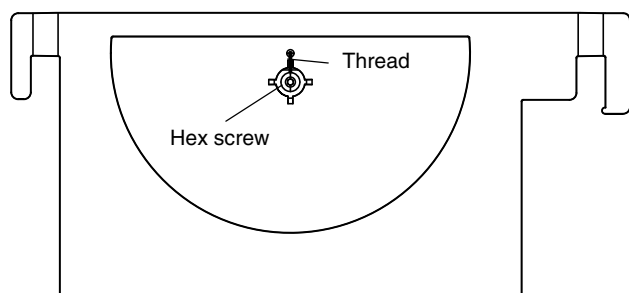

Loosen the hex screw in order to replace the thread. The table shows PASCO part numbers for replacement parts..

Table 1: PASCO Part Numbers

Part Number	Item
648-10296	Weight, brass
617-008	Thumbscrew
614-054	Square nut
610-270	Hex screw
699-011	Nylon thread

Limited Warranty

PASCO scientific warrants this product to be free from defects in materials and workmanship for a period of one year from the date of shipment to the customer. PASCO will repair or replace, at its option, any part of the product which is deemed to be defective in material or workmanship. This warranty does not cover damage to the product caused by abuse or improper use. Determination of whether a product failure is the result of a manufacturing defect or improper use by the customer shall be made solely by PASCO scientific. Responsibility for the return of equipment for warranty repair belongs to the customer. Equipment must be properly packed to prevent damage and shipped postage or freight prepaid. (Damage caused by improper packing of the equipment for return shipment will not be covered by the warranty.) Shipping costs for returning the equipment, after repair, will be paid by PASCO scientific.

Technical Support

For assistance with any PASCO product, contact PASCO at:

Address: PASCO scientific
 10101 Foothills Blvd.
 Roseville, CA 95747-7100

Phone: 916-786-3800 (worldwide)
 800-772-8700 (U.S.)

Fax: (916) 786-7565

Web: www.pasco.com

Email: support@pasco.com